

WarmStart® RTx Reverse Transcriptase



M0380S



50 reactions **15,000 U/ml** **Lot: 0011510**
RECOMBINANT **Store at -20°C** **Exp: 10/17**

Description: WarmStart RTx Reverse Transcriptase is a unique *in silico* designed RNA-directed DNA polymerase coupled with a reversibly-bound aptamer that inhibits RTx activity below 40°C. This enzyme can synthesize a complementary DNA strand initiating from a primer using RNA (cDNA synthesis) or single-stranded DNA as a template. RTx is a robust enzyme for RNA detection in amplification reactions and is particularly well suited for use in LAMP (loop-mediated isothermal amplification). The WarmStart property enables high throughput applications, room temperature setup, and increases the consistency and specificity of amplification reactions. RTx contains intact RNase H activity.

Source: An *E. coli* strain that carries the engineered RTx gene.

Application:

- RT-LAMP
- cDNA synthesis
- RT reactions requiring room temperature setup

Supplied in: 100 mM KCl, 10 mM Tris-HCl, 0.1 mM EDTA, 1 mM DTT, 50% glycerol, pH 7.4 @ 25°C

Reagents Supplied with Enzyme:

10X Isothermal Amplification Buffer, 100 mM MgSO₄

Reaction Conditions: 1X Isothermal Amplification Buffer, template, primer, dNTPs and 0.25–0.5 µl of WarmStart RTx Reverse Transcriptase in a reaction volume of 25 µl. Incubate at 50–55°C for cDNA synthesis or directly at 65°C for One-step RT-LAMP.

1X Isothermal Amplification Buffer:

20 mM Tris-HCl
10 mM (NH₄)₂SO₄
50 mM KCl
2 mM MgSO₄
0.1% Tween-20
pH 8.8 @ 25°C

Unit Definition: One unit is defined as the amount of enzyme that will incorporate 1 nmol of dTTP into acid-insoluble material in a total reaction volume of 50 µl in 20 minutes at 50°C using poly(rA)•oligo(dT)₁₈ as template.

Unit Assay Conditions: 20 mM Tris-HCl (pH 8.8), 10 mM (NH₄)₂SO₄, 50 mM KCl, 2 mM MgSO₄, 0.1% Tween-20, 0.5 mM dTTP, 0.4 mM poly(rA)•oligo(dT)₁₈.

Heat Inactivation: 10 minutes at 80°C

Typical RT-LAMP Protocol

COMPONENT	FINAL CONC./AMT.
10X Isothermal Amplification Buffer	1X
MgSO ₄ (100 mM)	6 mM (8 mM total)
dNTP Mix (10 mM)	1.4 mM
FIP/BIP Primers (25X)	1.6 µM
F3/B3 Primers (25X)	0.2 µM
LoopF/B Primers (25X)	0.4 µM
<i>Bst</i> 2.0 or <i>Bst</i> 2.0 WarmStart DNA Polymerase	8 units
WarmStart RTx Reverse Transcriptase	0.5 µl
Total Reaction Volume	25 µl

2. Mix by vortexing.
3. Place directly at 65°C.

Typical cDNA Synthesis Protocol:

COMPONENT	FINAL CONC./AMT.
10X Isothermal Amplification Buffer	1X
dNTP Mix (10 mM)	0.5 mM
Random Primer Mix	6 µM
RNase Inhibitor, Murine	20 units
WarmStart RTx Reverse Transcriptase	0.25 µl
Total Reaction Volume	20 µl

2. Mix by vortexing.
3. Incubate for 5 minutes at 25°C for annealing and 10 minutes at 55°C for synthesis.
4. Heat Inactivate WarmStart RTx Reverse Transcriptase by incubation at 80°C for 10 minutes.

Quality Control Assays

RNase Activity (Extended Digestion): A 10 µl reaction in 1X Isothermal Amplification Reaction Buffer containing 40 ng of labeled RNA and 15 units WarmStart RTx Reverse Transcriptase is incubated at 37°C. After incubation for 4 hours, > 90% of the substrate RNA remains intact as determined by agarose gel electrophoresis.

Endonuclease Activity (Nicking): A 50 µl reaction in 1X Isothermal Amplification Reaction Buffer containing 15 units WarmStart RTx Reverse Transcriptase with 1 µg ϕX174 RF I DNA for 4 hours at 37°C results in < 10% conversion to the RF II as determined by agarose gel electrophoresis.

Non-Specific DNase Activity (16-hour): A 50 µl reaction in 1X Isothermal Amplification Reaction Buffer containing 1 µg of Lambda DNA and 15 units of WarmStart RTx Reverse Transcriptase incubated for 16 hours at 37°C results in a DNA pattern free of detectable nuclease degradation as determined by agarose gel electrophoresis.

Exonuclease Activity (Radioactivity Release):

A 50 µl reaction in 1X Isothermal Amplification Reaction Buffer containing 1 µg of a mixture of single and double-stranded [³H] *E. coli* DNA and 15 units WarmStart RTx Reverse Transcriptase incubated for 4 hours at 37°C releases < 0.1% of the total radioactivity.

Protein Purity (SDS-PAGE): WarmStart RTx Reverse Transcriptase is ≥ 99% pure as determined by SDS-PAGE analysis using Coomassie Blue detection.

Companion Products Sold Separately:

Random Primer Mix
#S1330S 100 µl

RNase Inhibitor, Murine
#M0314S 3,000 units
#M0314L 15,000 units

Deoxynucleotide Solution Set
#N0446S 25 µmol each

Deoxynucleotide Solution Mix
#N0447S 8 µmol each
#N0447L 40 µmol each

Bst 2.0 DNA Polymerase
#M0537S 1,600 units
#M0537L 8,000 units

Bst 2.0 WarmStart DNA Polymerase
#M0538S 1,600 units
#M0538L 8,000 units

Isothermal Amplification Buffer Pack
#B0537S 6.0 ml



NEW ENGLAND BIOLABS®, THERMOPOL® and WARMSTART® are registered trademarks of New England Biolabs, Inc.

This product is intended for research purposes only. This product is not intended to be used for therapeutic or diagnostic purposes in humans or animals.

The purchase of NEB RTx products conveys to the purchaser the limited, nontransferable right to use the purchased products to perform reverse transcription loop-mediated isothermal amplification ("RT-LAMP") for research use only. LAMP is a patented technology belonging to Eiken Chemical Co., Ltd. and any use other than research may require a license from Eiken Chemical Co., Ltd. A patent is pending for NEB's RTx's product.